

GC/MS – Polymode Ionization – Pharmaceutical

## A Sample Prep-Free Analysis of Drug Molecules with “Smart IS+” and “SMCI+”

### Written by:

Chua Chun Kiang  
Tan Yu Jun  
Jackie Jackie  
Loo Lai Chin

### Abstract

The drug discovery and development journey is a very demanding process that would benefit from high throughput qualitative compound screening. As screening covers a wide range of chemical compounds, non-volatile compounds may be unsuitable for gas chromatography/mass spectrometry (GC/MS) analysis. This Technology Brief introduces a simple analysis technique that enables direct analysis of compounds with GC/MS. This is achieved by coupling a direct sample probe with innovative ion sources and reagent gas source capable of polymode ionization, which enable analysis of compounds in electron and chemical ionization analysis modes. The polymode ionization setup is expected to deliver speed and confidence in compound screening process with GC/MS.



*Keywords:*  
Polymode Ionization, Chemical Ionization,  
Drug Discovery, Drug Development,  
Direct Sampling

*Direct analysis of compounds such as inosine with GC-MS to accelerate drug discovery and development*

### Highlights

- “Smart IS+” and “SMCI+” setups direct qualitative GC/MS analysis of polar compounds such as drug molecules
- Choose between electron ionization and positive chemical ionization modes with “Smart IS+” setup or perform both ionization mode consecutively in a single GC/MS analysis
- Perform positive chemical ionization safely with “SMCI+” setup since it requires only common laboratory solvents as reagent gas

### Technologies Featured

#### Direct Sample Inlet



#### Smart EI/CI Ion Source (Smart IS)



#### Solvent Mediated Chemical Ionization (SMCI)



# 1. INTRODUCTION

The discovery and development of new drugs go through a series of lengthy processes that commonly begins in the synthetic laboratory. After the successful identification of a drug pharmacophore, the preparation of analogues demands a constant supply of starting materials. A rapid qualitative analysis method to confirm the successful synthesis of starting materials and thereafter the drug candidates would be important to increase the speed of drug discovery.

The usage of a direct probe as a sample inlet in GC/MS provides a rapid qualitative technique that eliminates the need for tedious sample preparation. It enables direct analysis of non-volatile, polar, and large molecules which typically requires multistep derivatization procedures for GC/MS analysis.

This article demonstrates the application of a direct probe in conjunction with a Smart EI/CI ion source (Smart IS) or solvent mediated chemical ionization (SMCI) unit for the analysis of drugs. The mass spectra of drug molecules generated by Smart IS and SMCI units will be evaluated.

# 2. EXPERIMENT

## 2.1 Analytical Setup

The analytical results in this report were generated using a Direct Sample Inlet (DI) probe in conjunction with a Smart IS or SMCI unit. The combination of DI with Smart IS or SMCI unit is hence known as “Smart IS+” and “SMCI+” in this article, respectively (Figure 1).

The DI probe is designed to be able to fit a miniature sample vial at its tip. The sample vial is thereafter placed close to the ion source and subsequently heated up according to a temperature program. The chemicals in the sample vial are hence volatilized and ionized in the ion source.

Smart IS is a 2-in-1 ion source that enables both electron ionization (EI) and positive chemical ionization (PCI) modes. PCI is achieved with the usage of isobutane gas as a reagent gas. Due to the simplicity of switching between two different ionization modes with Smart IS, the PCI mode attained with Smart IS is referred to as quick chemical ionization (QCI).

On the other hand, SMCI unit enables PCI mode with conventional PCI ion source and methanol as the reagent gas. Usage of methanol allows safe (i.e., it eliminates the use of flammable and toxic reagent gases such as methane, isobutane, and ammonia) and convenient adoption of PCI mode in routine GC/MS analysis.

A total of 7 drug molecules from two different classes, specifically, nucleic acid (i.e., adenosine, inosine) and antibiotic (i.e., sulfamethazine, sulfamethoxazole, furazolidone, furaltadone, tazobactam) were analyzed here.

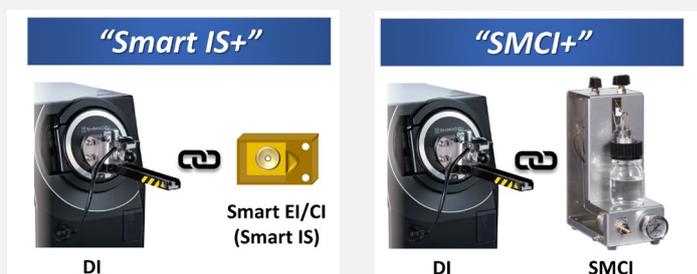


Figure 1. Polymode Ionization setup inclusive of “Smart IS+” and “SMCI+”

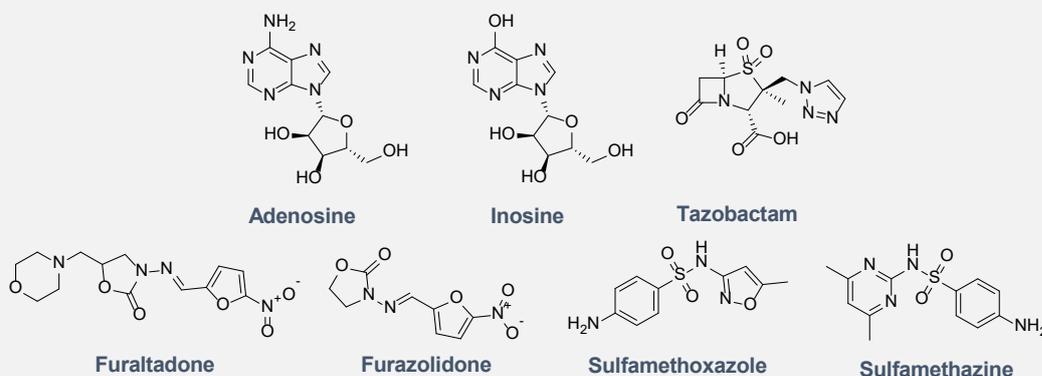


Figure 2. Drug molecules

## 2.2 Experimental Condition

Standard solutions of the drug molecules were prepared to a concentration of 5000 ppm in either water, acetone or DCM. 1  $\mu\text{L}$  of each standard solution was introduced into individual DI sample vial for analysis. The samples were left to dry before analysis.

The DI probe was heated at 20°C /min to 100°C, then 40°C /min to 450°C and held for 7 min. The ion source temperature was set to 230°C. Ionization mode used included EI, QCI (isobutane), and SMCI (methanol). Scan mode was performed in the range of  $m/z$  50-500 with a scan speed of 3333.

## 3. RESULTS AND DISCUSSION

### 3.1. “Mass Spectra of Nucleic Acids

Adenosine and inosine consist of a ribose bonded via a glycosidic bond. The presence of ribose rendered the molecules highly polar but was nevertheless analyzed directly with the DI probe. The EI mass spectra were successfully collected but the intensities of the molecular ion peaks were barely visible.


When analyzed with QCI mode, the distinctive  $[\text{M}+\text{H}]^+$  ion peak aid in the identification of the molecules. With SMCI mode, the  $[\text{M}+\text{H}]^+$  ion peak of adenosine was the base peak while that of inosine was distinctive but low in intensity.

### 3.2. Mass Spectra of Antibiotics

Sulfamethazine, sulfamethoxazole, furazolidone, and tazobactam are antibiotics for humans while furaltadone is a veterinary drug. The EI mass spectrum of furaltadone produced a clear fragment at  $m/z$  100 without any molecular ion signal, despite attaining a similarity index score of >80 against the NIST mass spectral library. With QCI and SMCI mode, obvious  $[\text{M}+\text{H}]^+$  ion peaks were observed, in which it appeared as the base peak in the latter.

Similarly, the molecular ion peak was not present in the EI mass spectrum of sulfamethazine, but the  $[\text{M}+\text{H}]^+$  ion peak was distinctive in both QCI and SMCI mass spectra. As for tazobactam, the EI mass spectrum did not consist of the molecular ion peak. In contrast, analysis with QCI mode resulted in a distinctive  $[\text{M}+\text{H}]^+$  ion peak at  $m/z$  301. The  $[\text{M}+\text{H}]^+$  ion peak was however not observed with SMCI mode.

While sulfamethoxazole and furazolidone produced EI mass spectra with obvious molecular ion peaks, both the QCI and SMCI mass spectra resulted in  $[\text{M}+\text{H}]^+$  ion peak as the base peak. With high similarity index scores for the EI mass spectra, the additional  $[\text{M}+\text{H}]^+$  ion peak information from the QCI and SMCI mass spectra further assisted in confirming their identities.

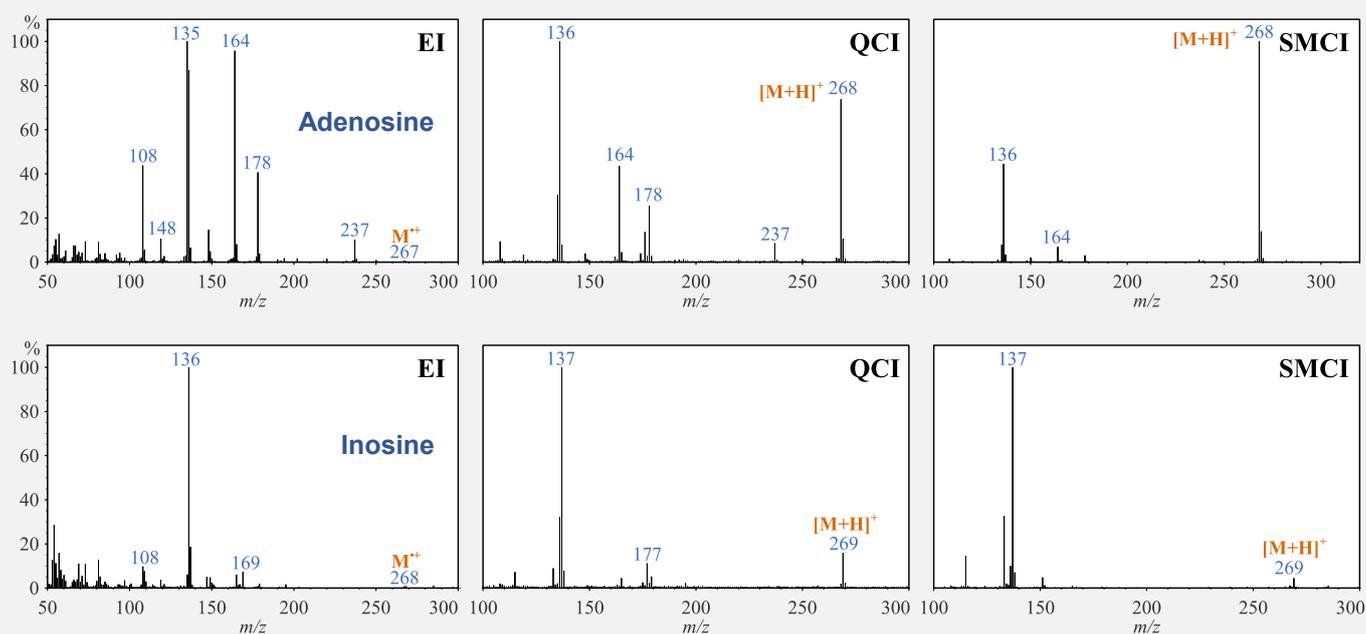

(continued on next page)

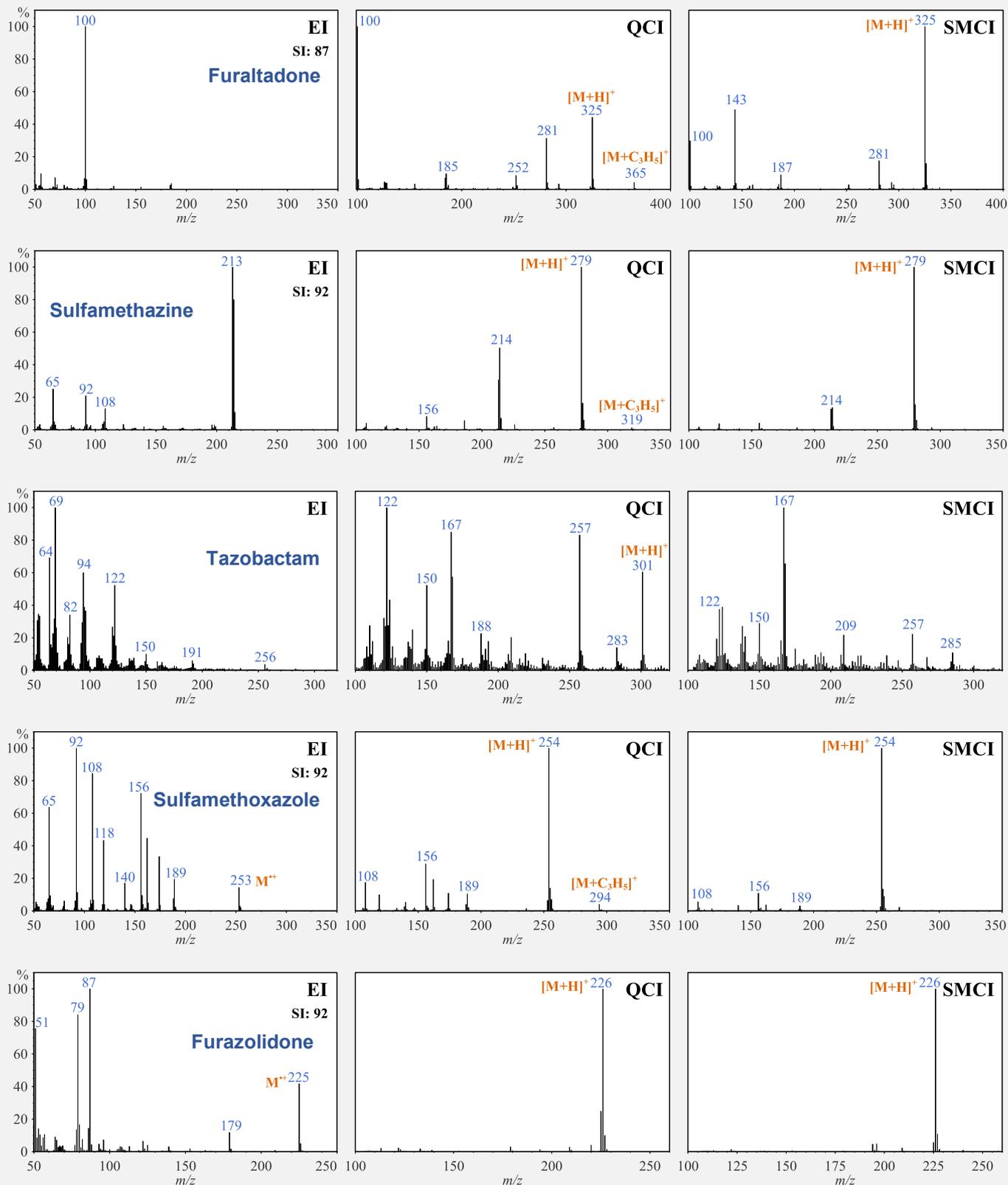


Figure 3. EI, QCI, and SMCI mass spectra of adenosine, inosine, furaltadone, sulfamethazine, tazobactam, sulfamethoxazole and furazolidone. SI: Library similarity index

## 4. CONCLUSION

The “Smart IS+” and newly introduced “SMCI+” enable a direct and quick qualitative analysis of drug molecules. Analysis of polar drug molecules with GC/MS can be accomplished directly and without the need for tedious derivatization steps.

The “Smart IS+” setup delivers convenience in switching between electron ionization and positive chemical ionization mode of analysis. In addition, a consecutive switch between EI and QCI mode can also be achieved within a single GC/MS run. On the other hand, the “SMCI+” setup delivers utmost convenience and safety to carry out positive chemical ionization since it utilizes methanol, which is a common laboratory solvent, as the reagent gas.

## READ MORE

### Direct Sample Inlet (DI)



The DI unit enables samples to be injected directly into the ion source without passing through the gas chromatograph. The DI probe can be heated up to 500°C under vacuum conditions. As a result, mass spectra of high polarity and high boiling point compounds can be generated easily. The analysis is performed by simply placing the sample in a sample vial that is secured in the DI probe.

[Learn More](#)

### Solvent Mediated Chemical Ionization (SMCI)

SMCI introduces headspace reagent gas from a bottle into the GC/MS ionization unit for use in positive or negative chemical ionization. Conventional chemical ionization techniques require the use of flammable reagent gas cylinders, but SMCI can be carried out with general organic solvents, such as methanol or acetonitrile, together with nitrogen or argon gas. This results in greater safety and lower running costs.



[Learn More](#)

### Ultra-High Sensitivity GC-MS



GCMS-TQ8050NX delivers the highest sensitivity for trace-level detection of genotoxic (mutagenic) impurities. Equipped with not only revolutionary OFF-AXIS ion optics for high ion transmission and outstanding noise elimination, but the GCMS-TQ8050NX also features a highly efficient detector capable of performing unprecedented quantitative analysis of ultra-trace amounts.

[Learn More](#)



Shimadzu (Asia Pacific) Pte. Ltd  
www.shimadzu.com.sg

**For Research Use Only. Not for use in diagnostic procedures.**

This publication may contain references to products that are not available in your country. Please contact us to check the availability of these products in your country.

The content of this publication shall not be reproduced, altered or sold for any commercial purpose without the written approval of Shimadzu.

Company names, products/service names and logos used in this publication are trademarks and trade names of Shimadzu Corporation, its subsidiaries or its affiliates, whether or not they are used with trademark symbol "TM" or "®".

Third-party trademarks and trade names may be used in this publication to refer to either the entities or their products/services, whether or not they are used with trademark symbol "TM" or "®".

Shimadzu disclaims any proprietary interest in trademarks and trade names other than its own.

The information contained herein is provided to you "as is" without warranty of any kind including without limitation warranties as to its accuracy or completeness. Shimadzu does not assume any responsibility or liability for any damage, whether direct or indirect, relating to the use of this publication. This publication is based upon the information available to Shimadzu on or before the date of publication, and subject to change without notice.